### **Nuance Software Release Notes**

Software Version: 3.0.2

This file provides information about Nuance software, version 3.0.2. Please read this information carefully as it supersedes all printed material or online help.

### **Overview:**

The Nuance software controls all Nuance / USB hardware.

## **System Requirements:**

- U.S. English Version of Microsoft<sup>©</sup> Windows XP (Service Pack 3 or higher) or U.S. English Version of Microsoft<sup>©</sup> Windows 7
- .NET 3.5 Service Pack 1
- Minimum monitor resolution: 1280x1024
- Minimum RAM: 2 GB

### Software Installation:

To install Nuance Software on Windows XP or Windows 7:

- 1. Log in as an administrator
- 2. Close all open software programs
- 3. Double click on Nuance's Setup.exe
- 4. Follow the installer's instructions

# Nuance Version 3.0.2

New Features:

- Install Nuance to PerkinElmer (ticket:873)
- Replace Caliper Life Sciences with PerkinElmer in the quick start guide (ticket:876)
- Replace Caliper Life Sciences with PerkinElmer in the manual (ticket:877)
- Update the splash screen (ticket:875)

Resolved Issues:

- Update the manual compute spectra algorithm (ticket:874)
- Have Nuance work with MetaMorph (ticket:878)

Known Issues:

• See 3.0.1.2 Know Issues

### Nuance Version 3.0.1.2

New Features:

None

Resolved Issues:

• The updated driver package was re-signed for Windows 7 64-bit compliance

Known Issues:

- Windows 7 only: Administrator privileges are required for running Nuance. When starting the software or opening files via Windows Explorer, Windows 7 will show a warning message stating that the Nuance software needs to access your hard drive. This is normal and will not cause problems.
- When connecting the hardware, administrator privileges are required to install the hardware drivers.

### **Documentation:**

• Complete documentation is accessible by selecting Help→Help within the Nuance software

- The User Manual is also available in .pdf format by selecting Help→View User Manual within the Nuance software or by accessing the Start menu (Caliper Life Sciences→Nuance→User Manual)
- The quick start guide is available in .pdf format by selecting Help→View Quick Start Guide within the Nuance software or by accessing the Start menu (Caliper Life Sciences→Nuance→Quick Start Guide)

## Support:

Contact Caliper Technical Support by phone, fax, or email from 8AM to 5PM Eastern Time.Phone:508.435.9761Toll Free (US only):877.522.2447Fax:508.435.3439Email:TechSupport@CaliperLS.comSupport Website:http://www.caliperls.com/support/contact-service-support.htm